

Lambda 365 Water Jacketed 8-Position Multi-Cell Holder Installation Instructions

This instruction sheet describes the installation of this accessory which is used with the Lambda 365 Spectrophotometer.

NOTE: *Read these instructions before you install this accessory.*

Contacting PerkinElmer

Supplies, replacement parts, and accessories can be ordered directly from PerkinElmer, using the part numbers.

See our website:

<http://perkinelmer.com>

PerkinElmer's catalog service offers a full selection of high-quality supplies.

To place an order for supplies and many replacement parts, request a free catalog, or ask for information:

If you are located within the U.S., call toll free 1-800-762-4000, 8 a.m. to 8 p.m. EST. Your order will be shipped promptly, usually within 24 hours.

If you are located outside of the U.S., call your local PerkinElmer sales or service office.

Features

- Suitable size for the standard cells
- Excellent durability
- Controls the temperature by circulation of liquid



Figure 1 Lambda 365 Water Jacketed 8-Position Multi-Cell Holder [P/N: N4101006]



PerkinElmer, 710 Bridgeport Avenue,
Shelton, CT06484-4794, U.S.A

Produced in the USA.

Dimensions and Specifications

Dimensions

Physical Characteristic		Specification	Comment
Outline	Height (mm)	143	
	Width (mm)	135	
	Depth (mm)	354	
Inner	Height (mm)	35	Suitable for the Standard Cell
	Width (mm)	14	
	Depth (mm)	12.6	
Weight (kg)		2.3	

Specifications

Physical Characteristic		Specification	Comment
Space between cells	(mm)	10.4	
Moving distance of one cell	(mm)	11.5	
Moving distance of eight cells	(mm)	80.5	
Moving time of one cell	(sec)	1.3	
Moving time of eight cells	(sec)	8.5	

Environmental Conditions for Operation

Specification	Comment
RT ~ 100 °C	Controls the temperature using Circulator

Configuration of Water Jacketed 8-Position Multi-Cell Holder

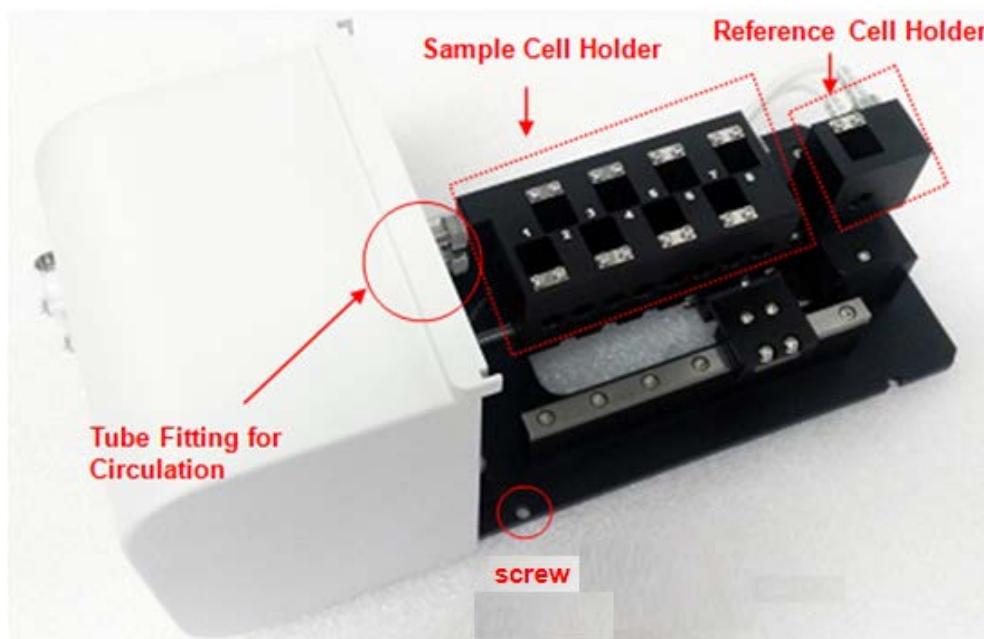


Figure 2 Water Jacketed 8-Position Multi-Cell Holder Configuration



- Used to fix a cell holder, a base plate or a front plate for Lambda 365.
- Spare screws (2ea) are enveloped with the accessory.

Figure 3 Phillips round headscrew with washer (M4 * 12L)



Figure 4. Male connector of quick coupling (2 each)

Installation

CAUTION *Make sure the instrument is turned off while installing this accessory.*
ATTENTION *Assurez-vous que l'instrument est éteint lors de l'installation de cet accessoire.*

1. Prepare the Lambda 365 Spectrophotometer to install this accessory.
2. Connect the power cord and the communication cable. **DO NOT** turn on the power of the instrument!
3. Remove the two Phillips round head screws with washer (M4*12L) to disassemble the existing cell holder and base plate



Figure 5. Location of the Phillips round head screws with washer

4. Pull out the cell holder and base plate by hand.



Figure 6. Pulling out the cell holder and base plate

5. Insert the Water Jacketed 8-Position Multi-Cell Holder in the sample compartment.



Figure 7. Inserting the Water Jacketed 8-Position Multi-Cell Holder into the sample compartment

6. Gently press the Water Jacketed 8-Position Multi-Cell Holder to connect the communication port (male) under the bottom of the Water Jacketed 8-Position Multi-Cell Holder to the port (female) in the sample compartment.

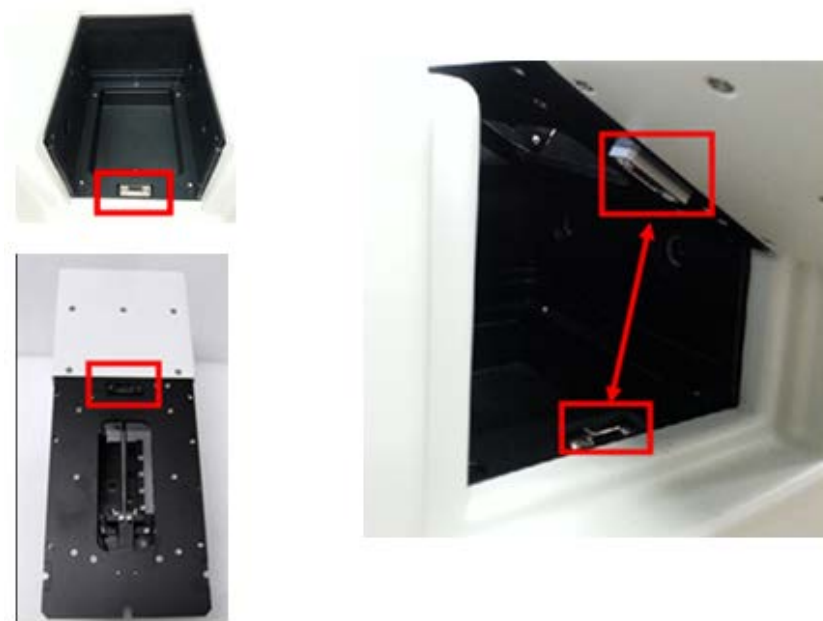


Figure 8. Connecting the communication ports

7. Tighten the Water Jacketed 8-Position Multi-Cell Holder in the sample compartment with the screws. (The red circles indicates the location of the screws in the figure).



Figure 9. Tighten the Water Jacketed 8-Position Multi-Cell Holder

8. Assemble the tube with the male connector of quick coupling.

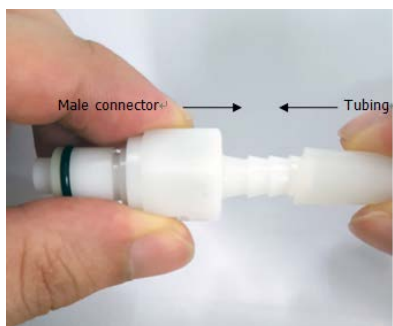


Figure 10. Assembling the tube with the male connector

NOTE: The tube size is recommended 6 mm-ID and 10 mm-OD silicon tube which has thermal resistance at least 100°C.

9. Connect the male connector with the tube to the quick coupling mount of the front plate.

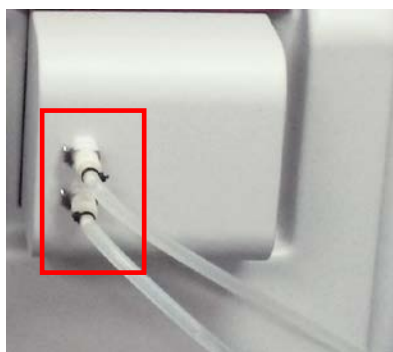


Figure 11. Connecting the male connector to the quick coupling

10. Turn on the power switch of the instruments.

Measurement


NOTE: Before executing System Self Test, the Water Jacketed 8-Position Multi-Cell Holder has to be installed first. Otherwise, the instrument can be damaged electrically and do not operate properly.

1. Install the Water Jacketed 8-Position Multi-Cell Holder.
2. Close the sample compartment cover and turn on the instrument.


CAUTION Do **not** start the UV Express software before finishing the initialization of grating.

NOTE: Start the System Self Test after warming up the system for at least 20 minutes.


3. Double-click on the **UV Express Software** folder and select one of the mode for starting.
4. The following message box will be shown. Empty the cell holder and close the lid firmly. Click **OK**.

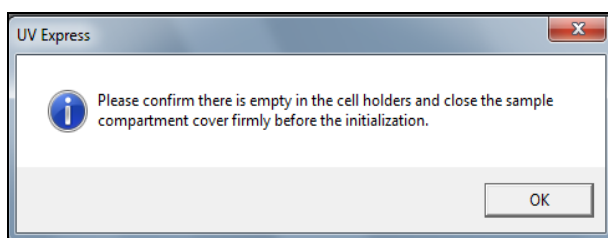


5. System Self Test starts. Click **OK** after finishing the **System Self Test**.
6. Select **New** to open a new window. Select **Experiment Type** and select **OK**.

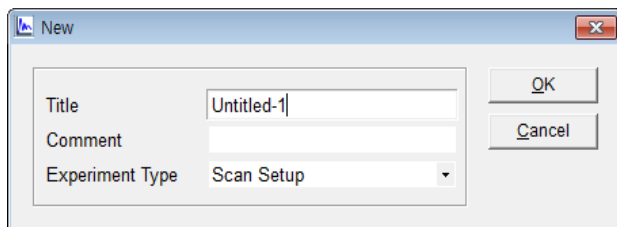



7. Open **Method** in the main menu.

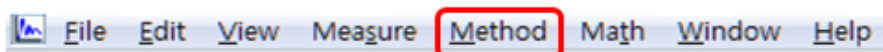

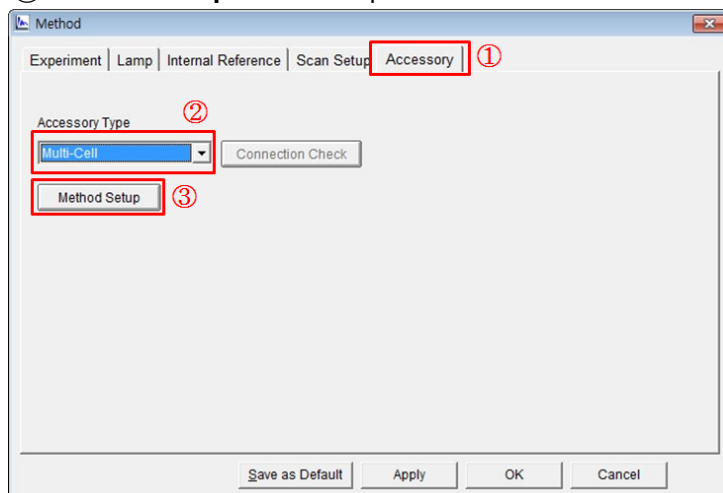


8. ① Select the **Accessory** tab and ② Select **Multi-cell** on **Accessory Type**, and then click the ③ **Method Setup** button to open the **Multicell** window.

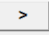

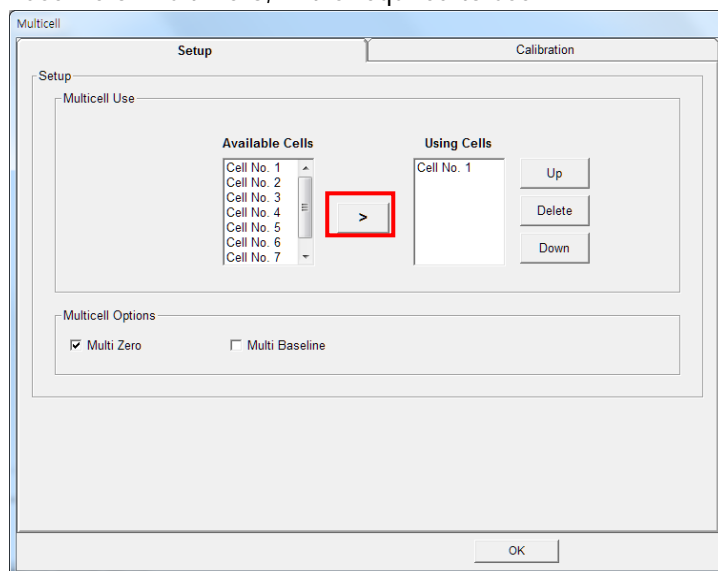9. Using the  button, select cells to be used among the available cells list. Check (✓) Multi Baseline or Multi Zero, if it is required to use.



- Available Cells:** Indicate cells that are available for measurement.
- Using Cells:** Shows the cell positions which are selected for measurement. Remove cells using  key and switch positions using  and  keys.
- Multicell Options**

Multi Zero: If checked (✓), zero will be measured all the selected positions. If not, zero will be only measured at the 1st cell position among the using cells.

Multi Baseline: If checked (✓), the baseline will be measured at each selected cell. If not checked, the baseline is only measured at the 1st cell position among the using cells.

NOTE: To prevent inadequate baseline (zero) measurement because of the difference in transmittance among cuvettes or any other reasons, Multi Baseline (Multi Zero) measurement is recommended.

10. Click **OK**.
11. After setting parameters for **Experiment, Lamp, Internal Reference and Scan Setup**, click **Apply** and then select **OK**.
12. After placing the reference solution into the selected Using Cell positions and the reference cell position of the Multicell holder, click the **Baseline** or **Zero** button to perform the Baseline or Zero function.


NOTE: *Baseline is required to run only one time when the first measurement is done, not every measurement. However, it should be done whenever the system is powered off/on, the wavelength is changed, SBW is changed or Reference sample is changed.*

13. Place the reference into the reference cell position and sample into the selected **Using cell** positions
14. Select the **Sample** button to measure.
15. After the measurement is finished, the results are displayed in the result window. Save or print the results as required.


Calibration of Multi-Cell Position

Calibrate the beam position of the Water Jacketed 8-Position Multi-Cell Holder when the Multicell is installed for the first time or beam position is incorrect.


1. Execute **Scan Setup** mode and ① click the **Accessory** tab of Method window. ② Select **Multi-cell** on **Accessory Type**, and then click the ③ **Method Setup** button to open the **Multicell** window.

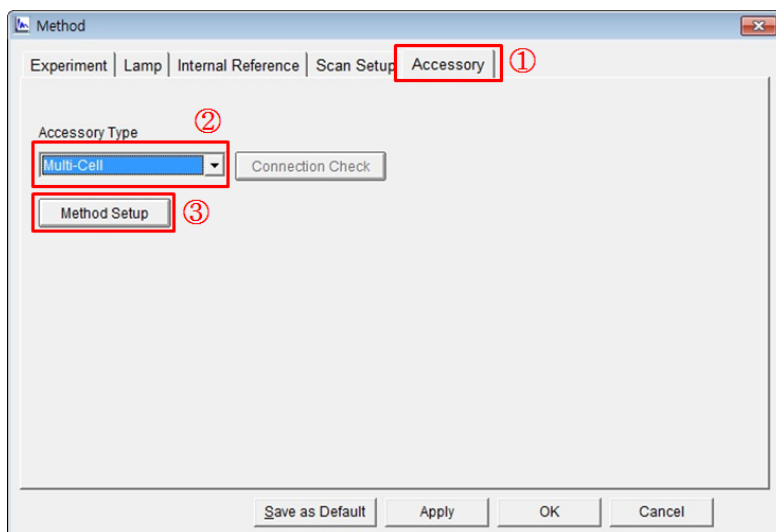


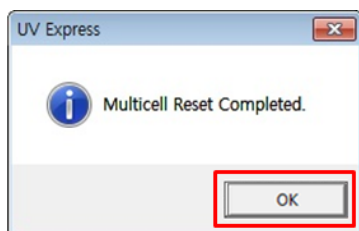
2. Select the **Calibration** tab in the **Multicell** window. The following dialog box will appear.

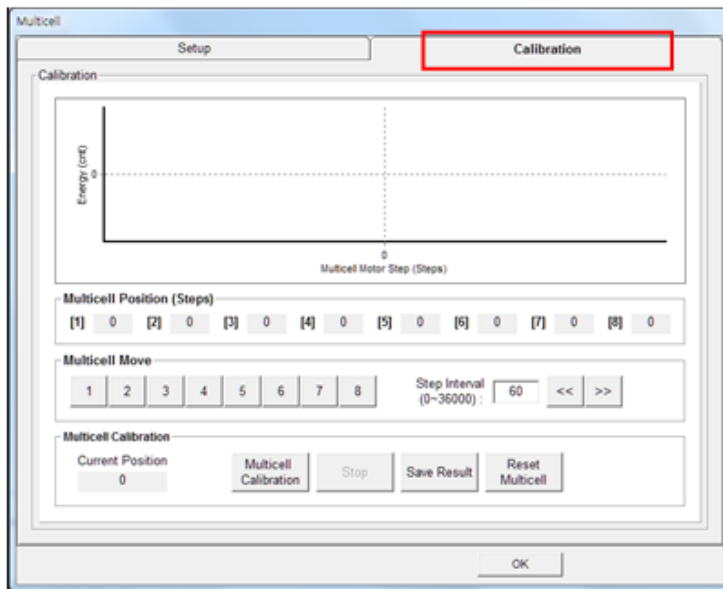



3. The functions of the Multicell Calibration are shown as follows.

Command	Function
MultiCell Position	Show saved steps about each cell position of the Multi-Cell.
Multicell Move	Used for moving Multi-Cell position as clicking buttons  Used for moving Multi-Cell position using << >> buttons by entered step.
Multicell Calibration	Used to perform the Multi-Cell calibration.
Stop	Used to stop the Multi-Cell calibration.
Save Result	Used to save the calibrated result.
Reset Multicell	Used to move to '0' step of Multi-Cell position.


4. Click on **Reset Multicell**  to format the Multi-Cell steps. Click **OK**.

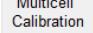


5. Select **MultiCell Calibration** , and then the following dialog box will appear. Remove all samples from the Multi-Cell holder (Empty the multicell holder).

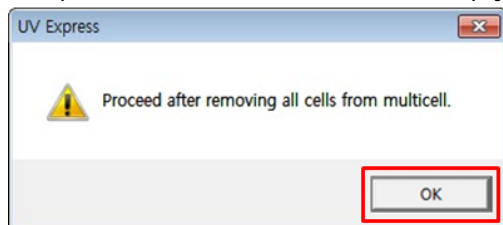



6. Select **OK**, and then the Multicell calibration will start. The current process of calibration will be shown in the main window.

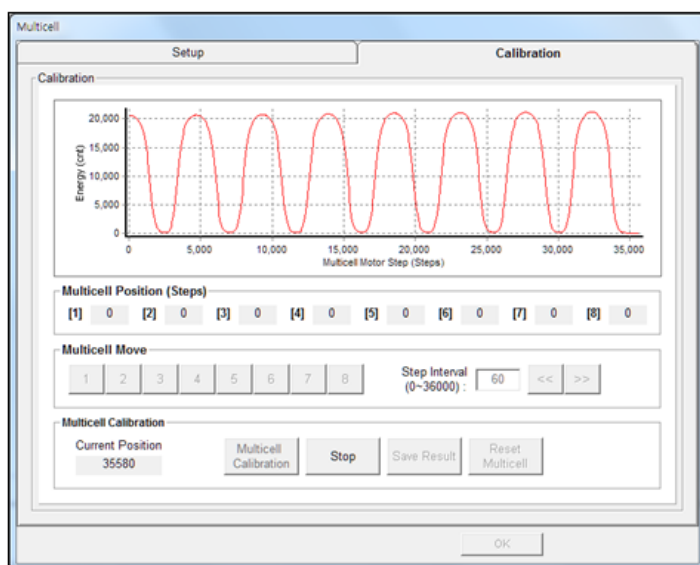

7. When calibration is finished, the following box appears. Select **OK**.

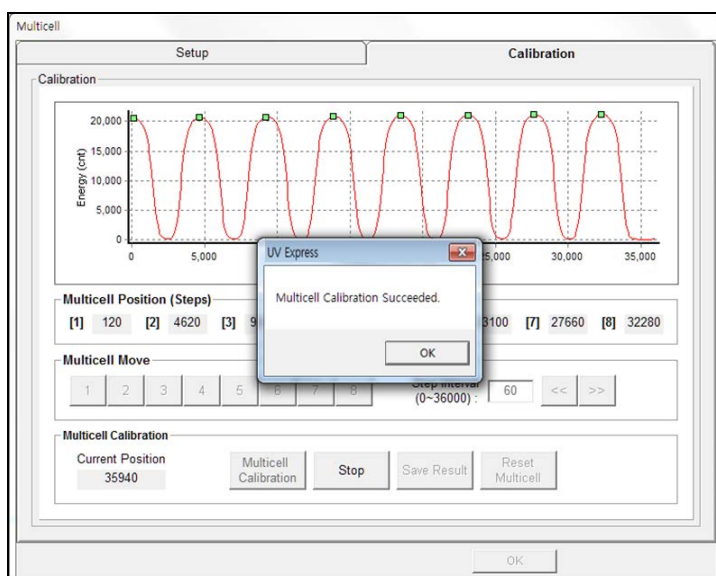

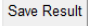8. Select **Save Result**  to save the result. If the following message box appears, select **OK**.

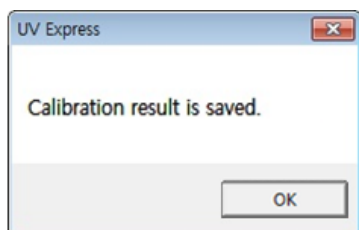

Troubleshooting

When liquid leakage occurs

1. When liquid is leaking at the fittings.
 - Using wrench, turn the fitting counter-clockwise to detach from the Water Jacketed Multi-Cell Holder.
 - Remove the Teflon which is wound around the fitting. Wind a length of new Teflon tape around the fitting.
 - Turn the fitting clockwise to attach it to the Water Jacketed Multi-Cell Holder.
2. When liquid is leaking between the quick coupling mount (or fitting) and the tube.
 - Replace the tube.
 - The replaced tube should be able to withstand at least 10 °C higher than the experimental temperature.
3. When liquid is leaking between the quick coupling mount and the Water Jacketed Multi-Cell Holder.
 - Contact PerkinElmer since the quick coupling mount needs to be replaced.

When constant temperature is not maintained

1. Check if the circulator is working normally.
 - Check if the temperature is maintained normally in the circulator.
 - Check if liquid is transferred normally to the Water Jacketed Multi-Cell Holder through the tube.
2. Check the tube connection.
 - Check if the tube is connected tightly.

When the tube frequently becomes detached

1. Replace the tube.
 - You should replace the tube as it is aged.

- The replaced tube should be able to withstand at least 10°C higher than the experimental temperature.
- 2. Fix with Teflon tape.
 - Firmly fix the tube by winding Teflon tape around the connection.

When the Multi-Cell does not move

1. Check with the connector port.
 - Check if the Multicell Holder connector is connected firmly to the Lambda 365.

When the intensity value is low

1. Recalibrate the Multi-Cell.
 - Recalibrate if the light beam does not go through the center of the cell holder.
2. Replace the lamp.
 - If the intensity value is still low after the recalibration, the lamps of the light source for measurement may be deteriorated and need replacing. Contact PerkinElmer for replacing the lamp.

When the Multi-Cell does not reset

If resetting fail, the photo interrupter switch inside the Multi-Cell will need replacing. Contact your PerkinElmer Service representative.