




ocm Anion Exchange Column				
	Mobile Phase Conditions			
Mobile Phase A	12mM NH4CO3			
Mobile Phase B	20mM (NH4)2HPO4			
HPLC Pump Conditions				
Separation Scheme	Gradient			
Pump Program	- Equal time: 7 min / A 100% - Run 1: 1 min / A 100% / Curve 0 - Run 2: 7 min / B 100%			
Pump Flow	2.0 mL/min			
Injection Volume	50 microliter			
ICP-MS Conditions				
Sample Introduction System	 Quartz Cyclonic Spray Chamber Glass Concentric Nebulizer 2.0 mm Quartz Injector			
Analyze Mode	AsO 91 / DRC Mode			
DRC Conditions	- O2 gas flow: 0.5ml/min - RPq: 0.45			
Dwell Time	500ms	_		

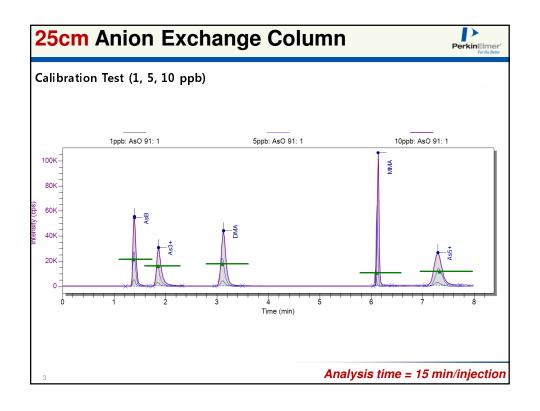

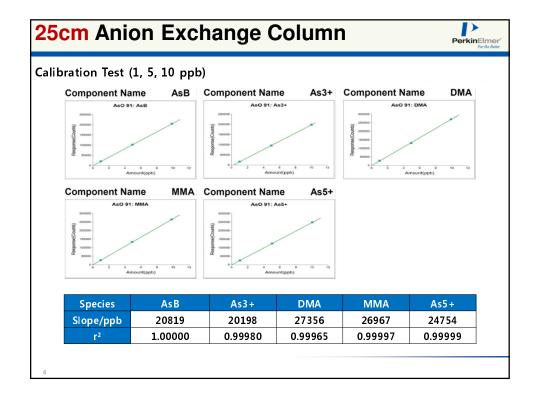


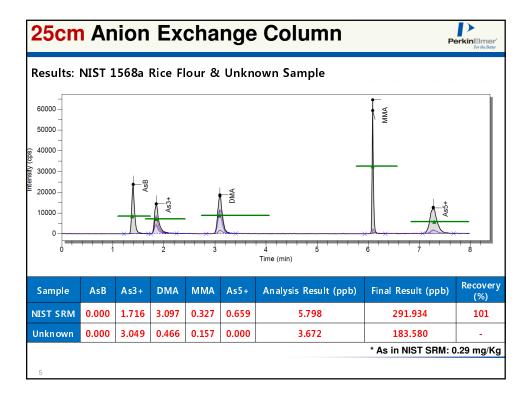

Conclusion



- 1. Very easy procedure of Mobile Phases
- 2. 15 Cm, 25 Cm Anion Exchange Column can be used for Speciation
 - 15 Cm Column: PKI Order Made
 - 25 Cm Column: Commercialized Products
- ➤ Disadvantage #1: Long analysis time (13~15min/injection)
- > Disadvantage #2: High base line = poorer Detection Limits

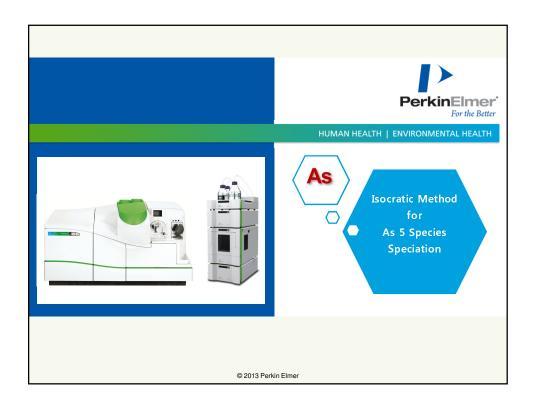

Preparing Your Lab.



- 1. Instruments: HPLC Speciation Kit, ICP-MS
- 2. Sample Preparation
 - Hot Block for 15 mL Tubes or Microwave, Centrifuge, Vortex Mixer
- 3. Bottles
 - 15 mL PE or PP Tube for Extraction / 2 mL PE or PP Tube for HPLC
 - 50 mL PP or Teflon Bottle for preparation for Organo-As Stock Standards
- 4. Column
 - Shiseido C18 MG II 5um, 250mm 4.6mm id
- 5. Stock Standards
 - Liquid Type As 3+ & 5+ (Purchased from ESI or SPEX or etc.)
 - Powder type "Arsenobetaine, Disodium methyl arsenate, Dimethylarsenic acid" (Purchased from Sigma Aldrich or etc.)

Preparing Your Lab. (Chemicals)



All chemicals purchased from Sigma-Aldrich

- 1.1-Octanesulfonic acid sodium salt (~98%) / C₈H₁₇NaO₃S / CAS: 6324-84-5
- 2.Malonic acid, ReagentPlus, 99% / C₃H₄O₄ / CAS: 141-82-2
- 3.Ammonium hydroxide, 5.0N in water / EC: 215-647-6
- 4. Analytical grade or Electronic grade Conc. HNO₃ (65%)
- ** Electronic grade Conc. HNO₃ from Dongwoo-Finechem in Korea
 - Density: 1.42 - %: 63~65%
 - Mole: 14.2 M
- ** 10% (v/v) HNO₃ & NH₄OH for pH adjustment of Mobile Phase

9

Sample Extraction Procedure: HPLC



- 1. Uncooked rice ground into a fine powder (like wheat flour)
- 2. Prepare the 0.03M (0.2% (v/v) HNO₃) solution → 0.1~1 Litter
- 3. 0.5 gram Rice Flour added to a 15mL Tube
- 4. 4.5 gram 0.2% HNO₃ solution added to a 15 mL Tube
- 5. Vortex (mixing) for 10 sec
- 6. Heating & Extraction with Hot Block @ 120 °C, 4 Hour
- 7. Cool down
- 8. Centrifuge @ 4,000 rpm, 30 min
- 9. Filter with PTFE Membrane (0.45 um)

10.0.1 mL Extracted Solution & 0.9 mL DIW added to a 2.0 mL HPLC Sample Tube

- 11. Vortex (mixing) for 10 sec
 - ** Microwave: Extract with Same condition # 6 using Extraction Mode in Microwave

Sample Digestion Procedure: Total As



- 1. Uncooked rice ground into a fine powder (like wheat flour)
- 2. 0.5 gram Rice Flour added to a 50 mL Tube
- 3. 4 mL Conc. HNO₃ & 1 mL Conc. H₂O₂ added to a 50 mL Tube
- 4. Vortex (mixing) for 10 sec
- 5. Heating with Hot Block @ 120 °C, 30 min
- 6. Cooling down
- 7. Bring to 50 mL with DIW

11

ICP-MS Method & STDs: Total As



- 1. Conditions for ICP-MS Method
 - Mass: AsO 91
 - DRC Gas: High purity Oxygen Gas
- CG / RPq: 0.5 / 0.65
- Internal Standard Element: Ir 193

2.STDs: 0.5, 1.0, 2.5 ppb (10 ppm ICP-MS Stock STD #3) 3.Internal STD: 5 ppb Ir (1,000 ppm Single Stock STD)

Preparation for Analysis (Stock STDs)



- 1. As 3+ & 5+ : Liquid Base
- 2. Organo-As: Powder Base / Need dilution with DIW
- 3. Preparation Method: Refer to below Equation

"As atom equivalent" = $\frac{\text{concentration of concentrated}}{\text{stock standard in mg/L}} \times \frac{\text{# As atoms per species molecule}}{\text{F.W. of As species}} \times 74.92 \text{ atomic wt. As}$

Arsenic Species	Formula Weight	Number of Arsenic Atoms per Molecule	Formula
Arsenobetaine	178.06	1	(CH ₃) ₃ AsCH ₂ COOH
Arsenocholine bromide	244.99	1	(CH ₃) ₃ AsCH ₂ CH ₂ OH·Br
Disodium methyl arsenate	291.9	1	CH ₃ AsO ₃ · 6H ₂ O
Trimethylarsine oxide	136.03	1	(CH ₃) ₃ AsO
Dimethylarsinic acid	138.01	1	(CH ₃) ₂ As(OH) ₂
Arsenic (V) oxide hydrate	229.84	2	As ₂ O ₅ · xH ₂ O
Arsenic (III) oxide	197.84	2	As ₂ O ₃

13

Preparation for Analysis (Stock STDs)



- 3. For example, prepare the 1,000 ppm As from AsB (50 mL)
 - 1,000 ppm = (? mg x 1 x 74.92) / 178.06 x 0.05 L

? = 118.833 mg (0.1188 g)

So, 0.1188 gram AsB added to a 50 mL Teflon Bottle, followed by DIW to 50 gram → Finally, 1,000 ppm As from AsB Stock Solution

- ** Other Organo-As: Same procedure with AsB
- 4. Mixed Stock Standard solution for Analysis: 10 ppm recommended

Dilute Solution: 0.2% HNO₃ Expiration Date: 1 month

5. Calibration Standard Solutions: 0.5, 2.5, 5.0 ppb recommended

Dilute Solution: DIW / Daily Preparation recommended


Preparation for Analysis




Prepare the Mobile Phase Solutions / Daily Preparation recommended


** pH & Mole Concentration are Very Important Spec


	HPLC Pump Conditions	
Separation Scheme	Isocratic	
Pump Program	- Equil time: 0.2 min / A 100% - Run: 3.5 min / A 100%	
Pump Flow	1.5 mL/min	
Injection Volume	50 microliter	
	ICP-MS Conditions	
Sample Introduction System	 Quartz Cyclonic Spray Chamber Glass Concentric Nebulizer 2.0 mm Quartz Injector 	
Analyze Mode	AsO 91 / DRC Mode	
DRC Conditions	- O_2 gas flow: 0.5ml/min - RPq: 0.45	
Dwell Time	500ms	

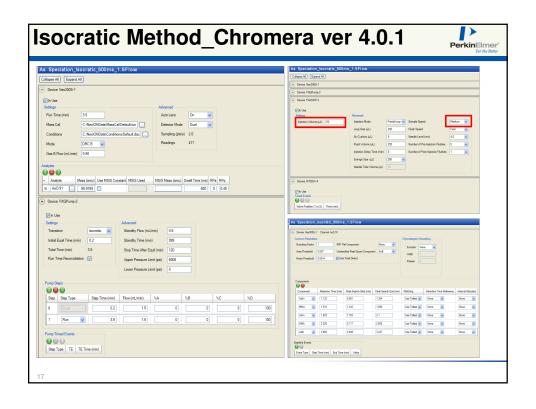


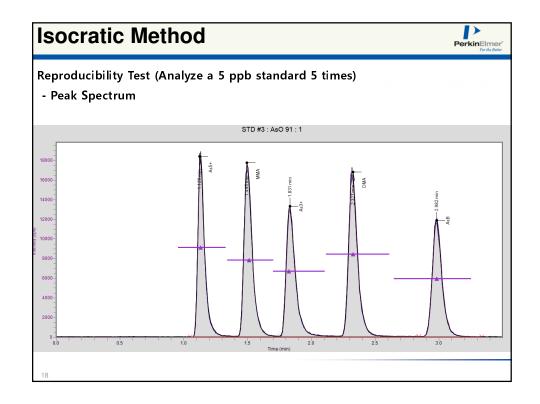


esults	=	est (An	alyze a	5 ppb	standa	ard 5 ti	mes)		
Species	RT #1	RT #:	2 RT	#3 R	tT #4	RT #5	AVG	SD	%RSD
As5+	1.126	1.120	5 1.1	34 1	134	1.134	1.131	0.004	0.387
ММА	1.495	1.49	5 1.5	04 1	.495	1.504	1.499	0.005	0.329
As3+	1.831	1.83	1.8	23 1	.831	1.831	1.829	0.004	0.196
DMA	2.318	2.32	7 2.3	18 2	2.327	2.318	2.322	0.005	0.212
AsB	2.982	2.982	2 2.9	74 2	2.982	2.982	2.980	0.004	0.120
Species	Results #1	Results #2	Results #3	Results #4	Results #	5 AVG	SD	%RSD	% Recovery
As5+	4.928	4.931	4.899	5.008	4.913	4.936	0.042	0.858	98.716
MMA	4.910	4.924	4.913	4.912	4.916	4.915	0.005	0.111	98.300
As3+	4.974	4.960	4.927	5.010	5.021	4.978	0.038	0.766	99.568
DMA	5.007	4.998	4.960	4.931	4.973	4.974	0.030	0.612	99.476
AsB	5.032	5.075	5.081	5.108	5.100	5.079	0.030	0.583	101.584

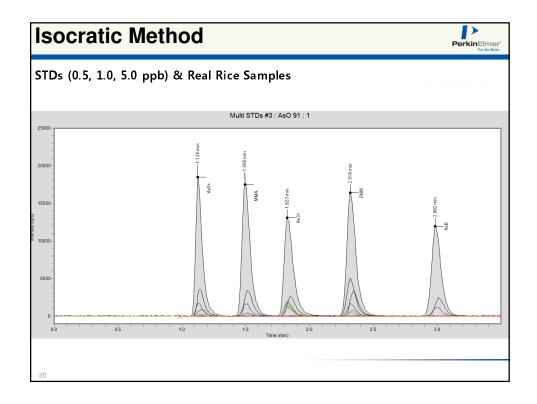

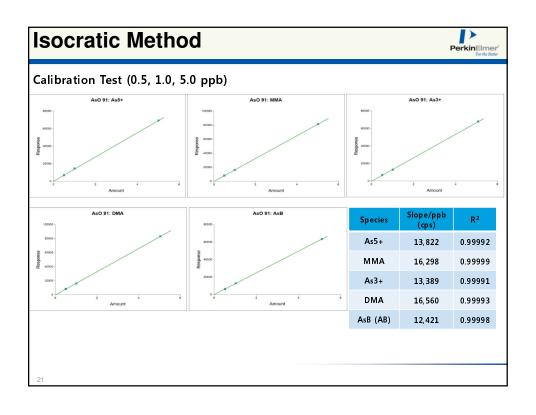

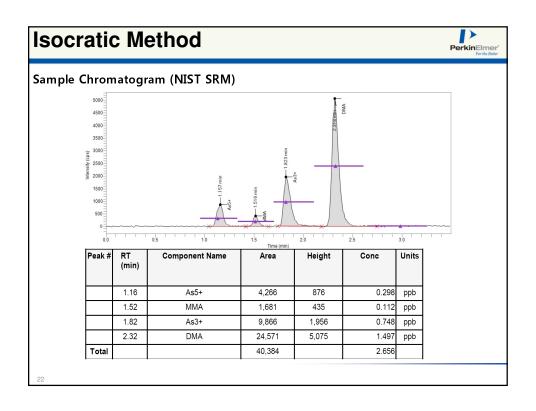

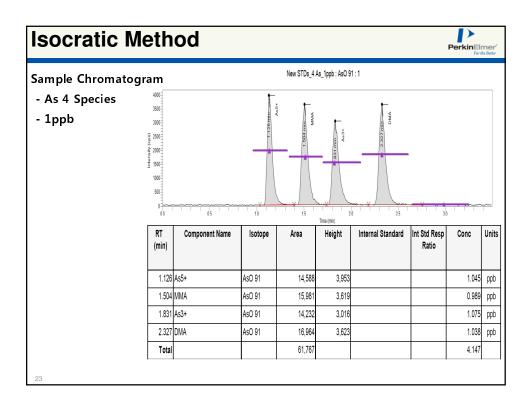

Final Results !!



Reference Material & Unknown Rice Sample

Results: ppb (ug/Kg)

#	Name -	Acid Digestion	Speciation	Speciation/Digestion	NIST SRM Recovery(%)	NIST SRM Recovery(%)
#	Name	AsO 91 (ppb)	AsO 91 (ppb)	Matching Rate(%)	Acid Digestion	Sp eciation
1	NIST SRM	287.36	265.60	92.43	99.09	91.59
2	Cooked Rice #1	68.03	73.40	107.89		
3	Cooked Rice #2	84.66	81.80	96.62		
6	Market Rice #1	106.11	116.80	110.07	NA	NA
5	Market Rice #2	95.46	91.00	95.33		
4	Market Rice #3	95.91	90.10	93.94		

- NIST SRM: 1568a Rice Flour (As in SRM: 290 ug/Kg)
- Cooked Rice: Steam Cooked Rice for Instant Food
- Market Rice #1: Cultivated from North Region of Korea
- Market Rice #2: Cultivated from Center Region of Korea
- Market Rice #3: Cultivated from South Region of Korea

Special Attention & Advantage




- 1. Accurate MP's pH & Mole Conc. are very important things
- 2. Final solution base of STDs for calibration is DIW
- 3. Base of Extraction Solution is 0.03M (0.2% (v/v) HNO₃) solution needed
- 4. Final Sample for Speciation is 10x dilution with DIW needed
- If used extraction solution's acidity is higher than recommended condition,
 all RT of peaks were delayed and increase the tailing → Poor Results
- 1. Advantage
 - Very Short Speciation Time: Totally 3.5 min
 - Very Low Baseline: AVG. 10~40 cps
 - Very Low Salt in Mobile Phase


PerkinEmer'
For the Better


HUMAN HEALTH | ENVIRONMENTAL HEALTH


Work on Animal feed


strument Pa	rameters			
	HPLC Pump Conditions			
Separation Scheme	Isocratic			
Pump Program	- Equil time: 0.5 min / A 100% - Run: 5 min / A 100%			
Column	Brownlee Validated C18 250mm 4.6mm 5 um N9303561			
Eluent	10mM Butanesulfonate 4mM Malonic + 4mM TMAH pH at 3 and 0.5% methanol			
Pump Flow	1.0 mL/min			
Injection Volume	50 microliter			
ICP-MS Conditions				
Sample Introduction System	 - Quartz Cyclonic Spray Chamber - Glass Concentric Nebulizer - 2.0 mm Quartz Injector 			
Analyze Mode	AsO 91 / DRC Mode			
DRC Conditions	- O ₂ gas flow: 0.8ml/min - RPq: 0.45			
Dwell Time	500ms			

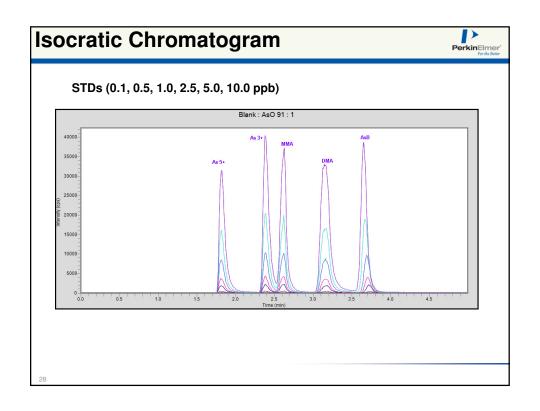


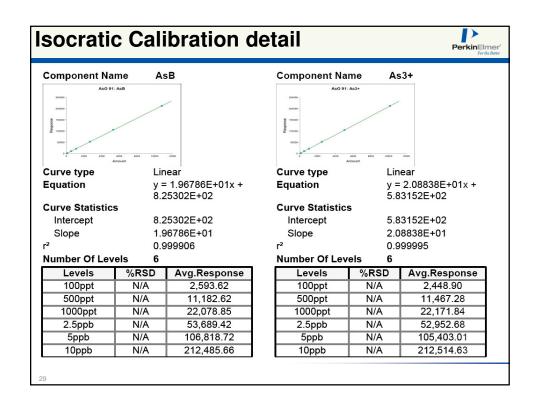


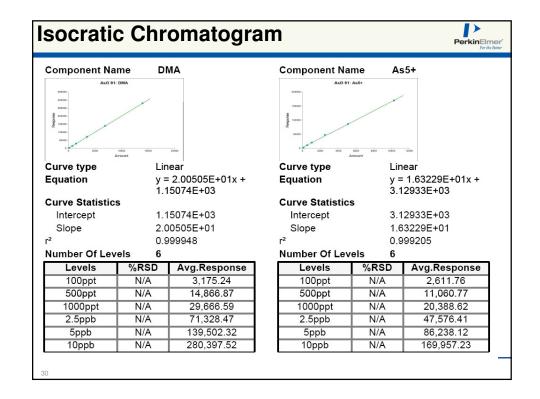

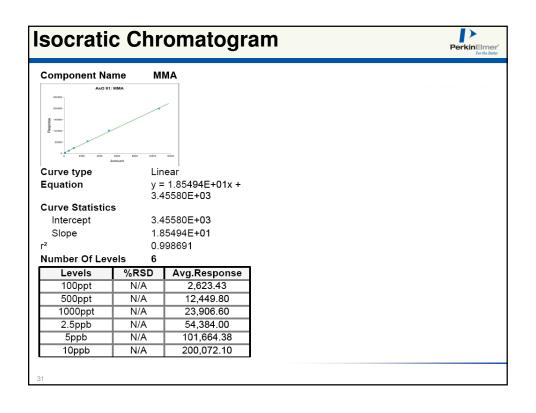


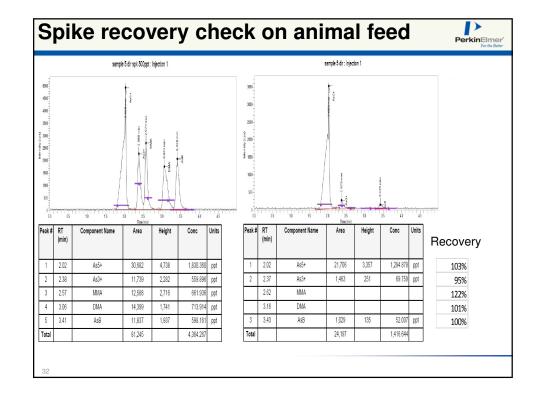

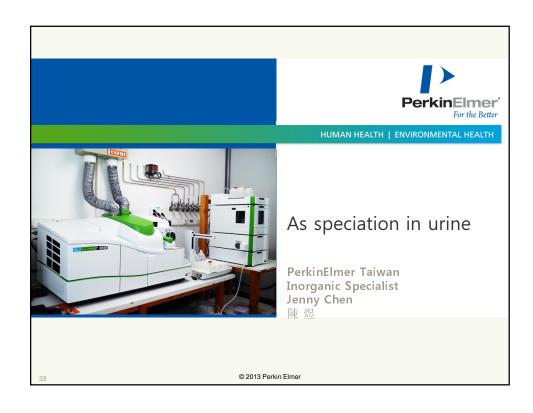

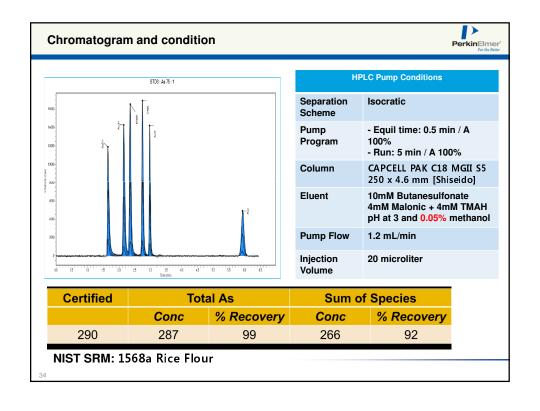

Calibration-AsIII and As Calibration curve: 0.1, 0.5, 1, 5, 10 ppb • Sample volume: 10 uL

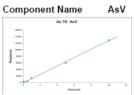


Curve type Linear Equation y = 1.28773E+03x -8.11066E+01

Curve Statistics

Intercept -8.11066E+01 Slope 1.28773E+03 0.997630

Number Of Levels

Levels	%RSD	Avg.Response		
STD1	N/A	201.24		
STD2	N/A	464.69		
STD3	4.237	1,360.68		
STD4	2.695	6,040.98		
STD5	1.872	12 940 17		

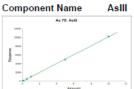

PerkinElme

 Curve type
 Linear

 Equation
 y = 1.01811E+03x +


Curve Statistics

Intercept 8.57067E-01 Slope 1.01811E+03 2 0.999010

Number Of Levels

Levels	%RSD	Avg.Response
STD1	N/A	142.70
STD2	4.401	522.70
STD3	5.263	1,090.15
STD4	0.897	4,912.55
STD5	1.541	10,263.49

8.57067E-01

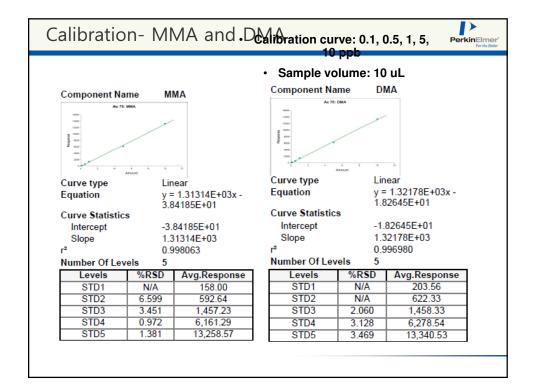

Calibration- AsB **Component Name** AsB Curve type Linear y = 1.10913E+03x -3.17514E+01 Equation Curve Statistics -3.17514E+01 Intercept 1.10913E+03 Slope 0.996850 **Number Of Levels** Avg.Response 135.74 Levels %RSD STD1 N/A 605.62 0.976 STD2 1,084.79 STD3 N/A STD4 1.368 5,277.43 STD5 4.400 11,173.01

