




AROMATICS HPLC

Environmental: BTEX Mix on Brownlee Analytical C18

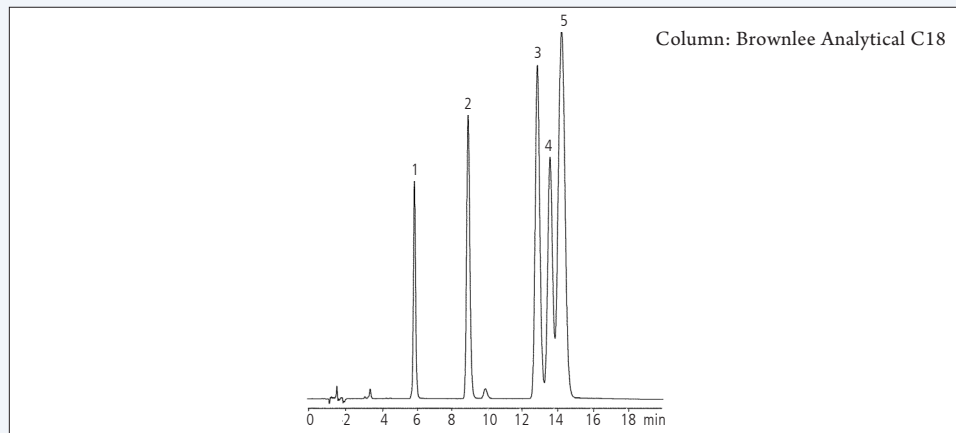

Peak List

1. benzene
2. toluene
3. o-xylene
4. ethylbenzene
5. m-xylene & p-xylene

Column

Column name: Brownlee Analytical C18
Part number: N9303513
Dimensions: 150 x 4.6 mm
Particle size: 5 μm
Pore size: 110Å

Conditions

Mobile phase: water:acetonitrile (50:50 v/v)
Flow: 1.0 mL/min
Temperature: ambient
Detector: UV @ 254 nm

Sample

Injection volume: 10 μL
Concentration: 200 $\mu\text{g/mL}$ each
Solvent: methanol

Notes

Recommended Guard Cartridge N9303677
Recommended Guard Holder N9303713

PerkinElmer, Inc.
940 Winter Street
Waltham, MA 02451 USA
P: (800) 762-4000 or
(+1) 203-925-4602
www.perkinelmer.com


PerkinElmer[®]
For the Better

For a complete listing of our global offices, visit www.perkinelmer.com/ContactUs

Copyright ©2013 PerkinElmer, Inc. All rights reserved. PerkinElmer[®] is a registered trademark of PerkinElmer, Inc. All other trademarks are the property of their respective owners.

007570A_06