




SUGARS/SWEETENERS

HPLC

Foods and Flavors: Sugar Analysis on Brownlee Polypore CA

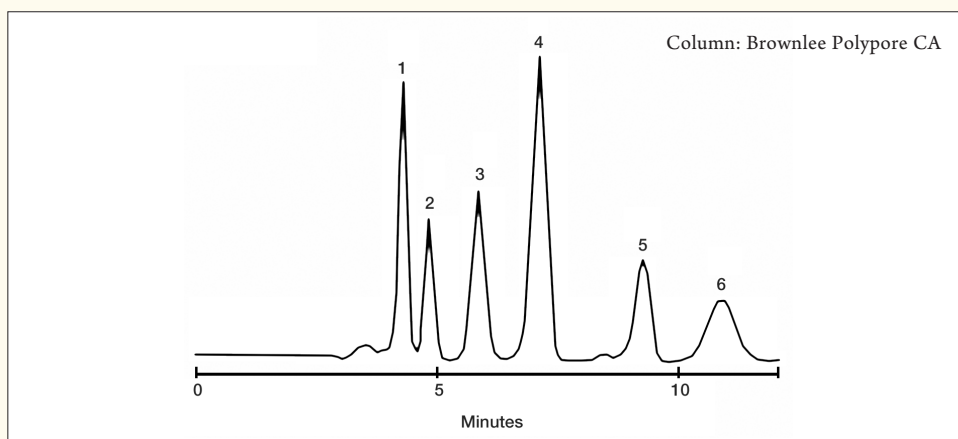

Peak List

1. maltotriose
2. maltose
3. glucose
4. fructose
5. mannitol
6. ribose

Column

Column name: Brownlee Polypore CA
Part number: 07110095
Dimensions: 220 x 4.6 mm i.d.
Particle Size: 10 μ m

Conditions

Mobile phase: water
Flow: 0.3 mL/min
Temperature: 85 °C
Detector: RI

Notes

Recommended Column Holder 07150015
Recommended Guard Cartridge 07110091
Recommended Guard Holder 07110013

PerkinElmer, Inc.
940 Winter Street
Waltham, MA 02451 USA
P: (800) 762-4000 or
(+1) 203-925-4602
www.perkinelmer.com

PerkinElmer
For the Better

For a complete listing of our global offices, visit www.perkinelmer.com/ContactUs

Copyright ©2013 PerkinElmer, Inc. All rights reserved. PerkinElmer® is a registered trademark of PerkinElmer, Inc. All other trademarks are the property of their respective owners.

007572A_37