

Lambda 465 Magnetic Film Holder Installation Instructions

This instruction sheet describes the installation of this accessory which is used with the Lambda 465 Spectrophotometer.

NOTE: *Read these instructions before you install this accessory.*

Contacting PerkinElmer

Supplies, replacement parts, and accessories can be ordered directly from PerkinElmer, using the part numbers.

See our website:

<http://perkinelmer.com>

PerkinElmer's catalog service offers a full selection of high-quality supplies.

To place an order for supplies and many replacement parts, request a free catalog, or ask for information:

If you are located within the U.S., call toll free 1-800-762-4000, 8 a.m. to 8 p.m. EST. Your order will be shipped promptly, usually within 24 hours.

If you are located outside of the U.S., call your local PerkinElmer sales or service office.

Features

- Easy to install
- Suitable size for the Standard Cell
- Easy to set Film
- Excellent durability



Figure 1 Lambda 465 Magnetic Film Holder [P/N: N4104007]

Dimensions and Specifications

Physical Characteristic		Specification	Comment
Outline	Height (mm)	81.5	
	Width (mm)	70	
	Depth (mm)	150	
Inner	Height (mm)	27.5	Suitable for the Standard Cell
	Width (mm)	12.5	
	Depth (mm)	12.5	
Weight (kg)		0.42	

Description

Configuration of the Magnetic Film Holder

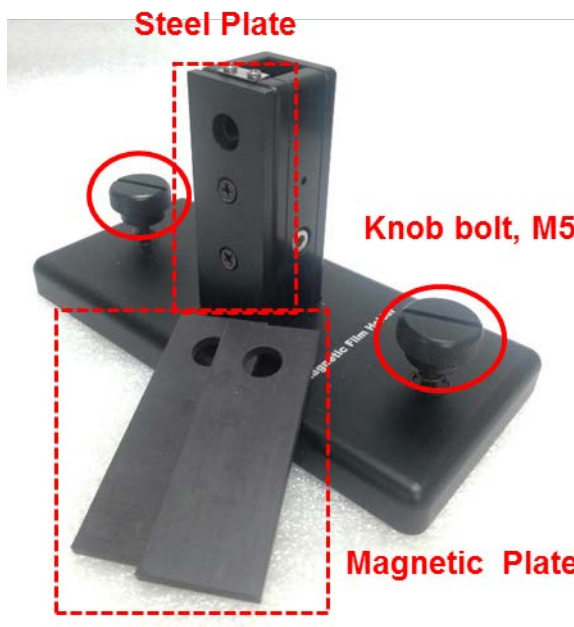


Figure 2 Lambda 465 Magnetic Film Holder

Installation

1. Disassemble the existing cell holder.

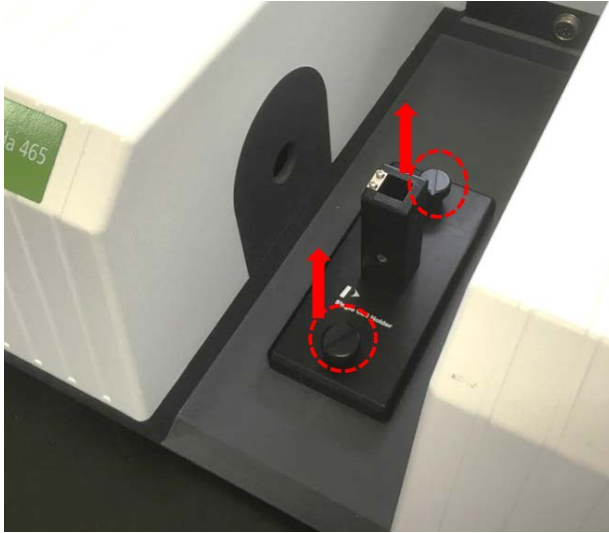


Figure 3 Location of the two knob bolts

2. Install the magnetic film holder by fastening the knob bolts.

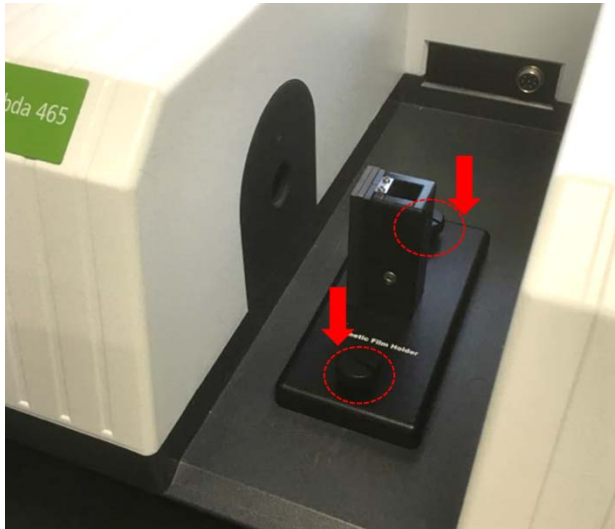


Figure 4 Installing the magnetic film holder

3. Connect the communication cable and the power cord, and then turn on the power of the Lambda 465.

